



Installation instructions of MODEM 1000 station



1. Installation

The station can be wall mounted (screws & plugs provided) or on a carriage, used horizontally or resting on a bench.



- Link the feeder pipe with the faucet by using the supplied adapter
- Remove the closing the tube output
- Adjust the flow rate to 60 liters/hour maximum

2. Change columns (K1000)

The saturation level of the station is visible: the saturated resin turns from GREEN to PURPLE.

- When the last column completely turned PURPLE, remove all columns from the support using the set of keys provided.
- Replace them by new columns.
- Replace the connections of input-output and those interconnecting columns.



3. Spare parts

DEM'O and MODEM 1000 replacement tip:

- elbow fitting : RC.DEMO

DEM'O and MODEM 1000 replacement pipe:

- polyethylene pipe : TP.DEMO

To prevent overpressure, close the blue water supply valve, upstream of the appliance, when not in use.